



Human Hepatic Stellate Cells (HHStC) Catalog #5300

Cell Specification

Hepatic stellate cells (HStC) are intralobular connective tissue cells presenting myofibroblast-like or lipocyte phenotypes. They participate in the homeostasis of liver extracellular matrix, repair, regeneration, fibrosis and control retinol metabolism, storage and release. Following liver injury, HStC transform into myofibroblast-like cells and are the major source of type I collagen in the fibrotic liver. Beyond these features, HStC have been implicated as regulators of hepatic microcirculation via cell contraction, and in disease states, in the pathogenesis of intrahepatic portal hypertension [1, 2]. Proliferation and migration of HStC and expression of chemokines are involved in the pathogenesis of liver inflammation and fibrogenesis. New insight into the molecular regulation of HStC activation will lead to therapeutic approaches in treatment of hepatic fibrosis, and could lead to reduced morbidity and mortality in patients with chronic liver injury.

HHStC from ScienCell Research Laboratories are isolated from human liver. HHStC are cryopreserved after purification and delivered frozen. Each vial contains $>5 \times 10^5$ cells in 1 ml volume. HHStC are characterized by immunofluorescence with antibodies specific to desmin and α -actin. HHStC are negative for HIV-1, HBV, HCV, mycoplasma, bacteria, yeast and fungi. HHStC are guaranteed to further expand for 15 population doublings under the conditions provided by ScienCell Research Laboratories.

Recommended Medium

It is recommended to use Stellate Cell Medium (SteCM, Cat. #5301) for the culturing of HHStC *in vitro*.

Product Use

HHStC are for research use only. They are not approved for human or animal use, or for application in *in vitro* diagnostic procedures.

Storage

Upon receiving, directly and immediately transfer the cells from dry ice to liquid nitrogen and keep the cells in liquid nitrogen until they are needed for experiments.

Shipping

Dry ice.

References

- [1] Reynaert H, Thompson MG, Thomas T, Geerts A. (2002) "Hepatic stellate cells: role in microcirculation and pathophysiology of portal hypertension." *Gut* 50:571-581.
- [2] Rockey DC. (2001) "Hepatic blood flow regulation by stellate cells in normal and injured liver." *Semin Liver Dis* 21(3):337-49.

Instructions for culturing cells

Caution: Cryopreserved cells are very delicate. Thaw the vial in a 37°C water bath and return the cells to culture as quickly as possible with minimal handling!

Initiating the culture:

1. Prepare a poly-L-lysine-coated culture vessel (2 $\mu\text{g}/\text{cm}^2$, T-75 flask is recommended). Add 10 ml of sterile water to a T-75 flask and then add 15 μl of poly-L-lysine stock solution (10 mg/ml, Cat. #0413). Leave the vessel in a 37°C incubator overnight (or for a minimum of one hour).
2. Prepare complete medium. Decontaminate the external surfaces of medium bottle and medium supplement tubes with 70% ethanol and transfer them to a sterile field. Aseptically transfer supplement to the basal medium with a pipette. Rinse the supplement tube with medium to recover the entire volume.
3. Rinse the poly-L-lysine-coated vessel twice with sterile water and then add 15 ml of complete medium. Leave the vessel in the sterile field and proceed to thaw the cryopreserved cells.
4. Place the frozen vial in a 37°C water bath. Hold and rotate the vial gently until the contents completely thaw. Promptly remove the vial from the water bath, wipe it down with 70% ethanol, and transfer it to the sterile field.
5. Carefully remove the cap without touching the interior threads. Gently resuspend and dispense the contents of the vial into the equilibrated, poly-L-lysine-coated culture vessel. A seeding density of 5,000 cells/ cm^2 is recommended.

Note: Dilution and centrifugation of cells after thawing are not recommended since these actions are more harmful to the cells than the effect of residual DMSO in the culture. It is also important that cells are plated in poly-L-lysine-coated culture vessels to promote cell attachment.

6. Replace the cap or lid of the culture vessel and gently rock the vessel to distribute the cells evenly. Loosen cap, if necessary, to allow gas exchange.
7. Return the culture vessel to the incubator.
8. For best results, do not disturb the culture for at least 16 hours after the culture has been initiated. Refresh culture medium the next day to remove residual DMSO and unattached cells, then every other day thereafter.

Maintaining the culture:

1. Refresh supplemented culture medium the next morning after establishing a culture from cryopreserved cells.
2. Change the medium every three days thereafter, until the culture is approximately 70% confluent.

3. Once the culture reaches 70% confluency, change medium every other day until the culture is approximately 90% confluent.

Subculturing:

1. Subculture when the culture reaches 90% confluency or above.
2. Prepare poly-L-lysine-coated culture vessels ($2 \mu\text{g}/\text{cm}^2$) one day before subculture.
3. Warm complete medium, trypsin/EDTA solution (T/E, Cat. #0103), T/E neutralization solution (TNS, Cat. #0113), and DPBS (Ca^{++} - and Mg^{++} -free, Cat. #0303) to **room temperature**. We do not recommend warming reagents and medium in a 37°C water bath prior to use.
4. Rinse the cells with DPBS.
5. Add 8 ml of DPBS and then 2 ml of T/E solution into flask (in the case of a T-75 flask). Gently rock the flask to ensure complete coverage of cells by T/E solution. Incubate the flask in a 37°C incubator for 1 to 2 minutes or until cells completely round up. Use a microscope to monitor the change in cell morphology.
6. During incubation, prepare a 50 ml conical centrifuge tube with 5 ml of fetal bovine serum (FBS, Cat. #0500).
7. Transfer T/E solution from the flask to the 50 ml centrifuge tube (a small percent of cells may detach) and continue to incubate the flask at 37°C for another 1 to 2 minutes (no solution in the flask at this moment).
8. At the end of incubation, gently tap the side of the flask to dislodge cells from the surface. Check under a microscope to make sure that all cells detach.
9. Add 5 ml of TNS solution to the flask and transfer detached cells to the 50 ml centrifuge tube. Rinse the flask with another 5 ml of TNS to collect the residual cells.
10. Examine the flask under a microscope for a successful cell harvest by looking at the number of cells being left behind; there should be less than 5%.
Note: Use ScienCell T/E solution that is optimized to minimize cell damages due to over trypsinization.
11. Centrifuge the 50 ml centrifuge tube at 1000 rpm for 5 minutes. Resuspend cells in culture medium.
12. Count and plate cells in a new poly-L-lysine-coated culture vessel with the recommended cell density.

Caution: Handling human derived products is potentially biohazardous. Although each cell strain tests negative for HIV, HBV and HCV DNA, diagnostic tests are not necessarily 100% accurate, therefore, proper precautions must be taken to avoid inadvertent exposure. Always wear gloves and safety glasses when working with these materials. Never mouth pipette. We recommend following the universal procedures for handling products of human origin as the minimum precaution against contamination [1].

[1] Grizzle WE, Polt S. (1988) "Guidelines to avoid personal contamination by infective agents in research laboratories that use human tissues." *J Tissue Culture Methods*. 11: 191-9.